



# Master Touch series water purification system

Master Touch series is optimized and upgraded on the basis of Master series, which is the sole leading brand of Chinese Good Instrument in lab water area. It is the most representative products in Chinese lab water market.

## ◆ Features And Advantages

5.0 inch colorful high-resolution touch screen (16:9) controlling system, achieve finger-touch new experience

3 way online water quality sensor, detect the quality of feed water, RO water, deionized water, or ultrapure water respectively. And warn once water quality's standard exceeding

2 way flow sensor, achieve quantified dispensing of RO water, deionized water, or ultrapure water

System sanitizing procedure, achieve the disinfection of ultrapure water's tube and valve

System circulation function, achieve ultrapure water's circulation to keep top quality of ultrapure water

All Cartridges replacing alarm function, based on time, or water quality, show cartridges' used and residual life

Multiple alarm function: no feed water, full water, water quality's standard exceeding, and cartridge life ending

Auto self-flushing of RO membrane function (interval and continuous time setting), extend RO membrane's life

Auto running data storing function with built-in SD card, and data can be exported through the USB interface

Comprehensive Information query and management function, master system status, water quality, cartridges usage and alarm information

System time setting (year/month/day/hour/minute), timing standby (0-60 minute), and timing shutdown (0-24 hour) function

Level II password, protect all the parameters setting, and prohibit any unauthorized setting change

2 kind of pure water tank (liquid level PE tank and pressure tank). Also external tanks is optional

Whole plastic shell with high-strength, avoid rusting and keep clean, to meet GLP standard

3 door and easy-to-replacing cartridge design, convenient to maintain system and replace cartridges

Tube and adapter with NSF authorization and top quality, reduce TOC level and assure ultrapure water's quality

Optimized pretreatment (including PP fiber, KDF and active carbon cartridge), effectively protect RO membrane

RO module with DOW's membrane, ensure long life, stable operation and high desalinization rate

4 in 1 ultrapure cartridge (can be divided to 4 independent cartridge) with DOW's top polishing resin, ensure ultrapure water's quality up to 18.2 MΩ.cm, with the lowest TOC level

Double wavelength (185&254nm) ultraviolet lamp module, restrain bacteria's increase and reduce TOC


MWCO 5000D ultrafiltration module, effectively eliminate endotoxin, and suitable for precise cell cultivating and IVF

(0.45±0.1)μm double layer PES terminal disinfection filter, assure that terminal pure water is absolutely axenic



# Master Touch-Q series deionized water system (Tap water inlet)

## ◆ Flow Schematic

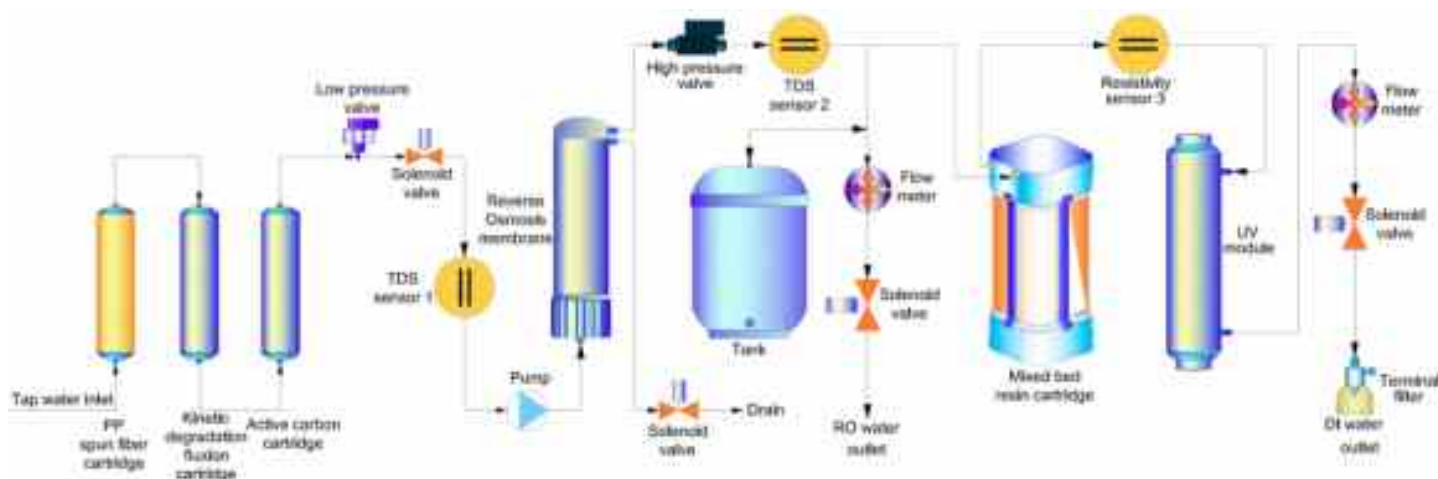

## ◆ Specifications

Model	Master Touch-Q15	Master Touch-Q15UT	Master Touch-Q30	Master Touch-Q30UT
Output(25°C)*	15 liters/hour		30 liters/hour	
Flow rate	Up to 2 liters/minute (with pressure tank)			
Pure water outlet	2: reverse osmosis water, deionized water			
Deionized water quality				
Resistivity	16-18.2MΩ.cm			
Bacteria	N/A	<0.1cfu/ml	N/A	<0.1cfu/ml
Particle(>0.1µm)	N/A	<1/ml	N/A	<1/ml
RO water quality				
Ion rejection rate	96%-99% (new RO membrane)			
Organic rejection rate	>99%, when MW>200 Dalton			
Particles and bacteria rejection rate	>99%			
Feed water requirements	Tap water, temperature:5-45°C,pressure:1.0-4.0Kgf/cm <sup>2</sup>			
Dimension and weight	Length×Width×Height:500×360×540mm / Weight: about 20Kg			
Electrical requirements	AC110-240V, 50/60Hz			
Power	120W			
Standard configuration	Main body (Including 1 set of cartridge)+15 liters pressure tank			

### Remarks:

\*The value will be influenced by temperature and feed water's quality.





Particles and bacteria rejection rate	>99%
Feed water requirements	Tap water, temperature:5-45°C,pressure:1.0-4.0Kgf/cm <sup>2</sup>
Dimension and weight	Length×Width×Height:500×360×540mm / Weight: about 20Kg
Electrical requirements	AC110-240V, 50/60Hz
Power	120W
Standard configuration	Main body (Including 1 set of cartridge)+15 liters pressure tank

**Remarks:**

\*The value will be influenced by temperature and feed water's quality.






# Master Touch-D series ultrapure water system (Distilled water inlet)

- With 5.0 inch touch screen system, 3 way water quality sensor, 2 way flow sensor for quantified dispensing and 1 pump, Master Touch-D series ultrapure water system is superior choice of ultrapure water for high grade experiments.
- With pure water or distilled water inlet, its output is up to 2 liters/minute. It can produce deionized water and ultrapure water. The deionized water's resistivity is above 5MΩ.cm, and the ultrapure water's resistivity absolutely reaches to 18.2MΩ.cm. It completely meets the highest grade I standard of ASTM, CAP, CLSI, EP and USP.

## ◆ Flow Schematic

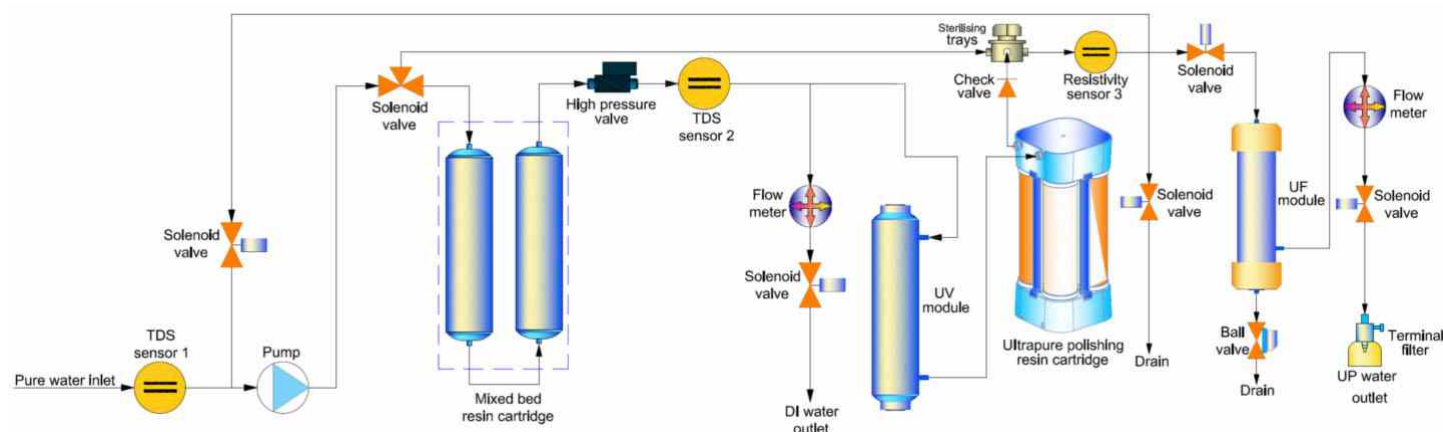

## ◆ Specifications

Model	Standard	Eliminating endotoxin	Low TOC	Synthesizing
	Master Touch-D	Master Touch-DUF	Master Touch-DUV	Master Touch-DUVF
Output	Up to 2 liters/minute (less output with UF cartridge)			
Pure water outlet	2: deionized water, ultrapure water			
Ultrapure water quality				
Resistivity(25°C)	18.2MΩ.cm			
TOC*	<10ppb	<10ppb	<3ppb	<3ppb
Bacteria	<0.1cfu/ml			
Particle(>0.1μm)	<1/ml			
Endotoxin	N/A	<0.001Eu/ml	N/A	<0.001Eu/ml
RNases	N/A	<0.01ng/ml	N/A	<0.01ng/ml
DNases	N/A	<4pg/μl	N/A	<4pg/μl



Deionized water quality	
Resistivity(25°C)	>5MΩ.cm
Feed water requirements	RO water, distilled water, deionized water, 5-45°C, 1atm*
Dimension and weight	Length×Width×Height:500×360×540mm / Weight: about 18Kg
Electrical requirements	AC110-240V, 50/60Hz
Power	120W
Standard configuration	Main body (Including 1 set of cartridge)

**Remarks:**

\*The value will be influenced by temperature and feed water's quality.

